



Instructions for using Powder Thief

Thoroughly read this sheet before using the device

General:

The Powder Thief is designed for use with free-flowing powders and granules.

Assembly:

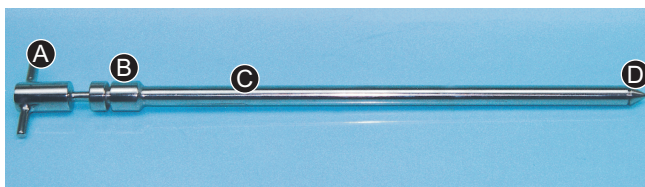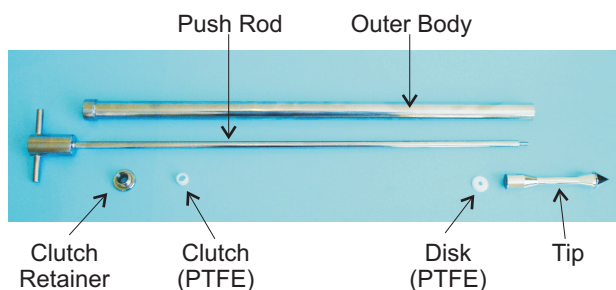

- (A) Handle
- (B) Clutch Housing
- (C) Body
- (D) Detachable Tip




Detachable Tips

1. Push the Clutch Retainer and then the Clutch up the push rod. Ensure that the clutch is correctly orientated




2. Insert the Push Rod into the Outer Body - lightly screw in the Clutch Retainer
3. Push the Disk over the threaded end of the Push Rod

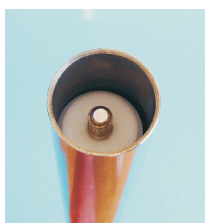

4. Screw the Tip into position - ensure that it is screwed in as far as it will go (about 6 turns)
5. Adjust the Clutch Retainer to give sufficient drag to hold the push rod in place when the sampler is held vertically

Operation:

1. Insert the sampler into the product, ensure that the tip is inside the sampler body.
2. At the required depth pull up the body to expose the tip. Powder will flow in around the tip.
3. Push down body of the sampler to trap the sample.
4. Withdraw sampler.
5. Pull up body to release the sample.

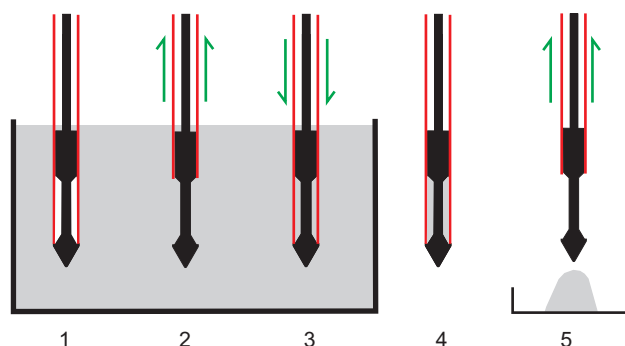

Cleaning:

The sampler should be stripped down to its component parts and then cleaned according to the usual cleaning methods that are employed for items such as general dispensing ware, tablet punches etc.

Ensure that the all component parts are thoroughly dried after washing.

Hand Wash ✓	Washing Machine ✓	Ultra-sonic bath ✓	Autoclave 121°C ✓
-------------	-------------------	--------------------	-------------------

IMPORTANT: ENSURE SAMPLER IS CLEANED THOROUGHLY BEFORE INITIAL USE

Whilst every care has been made during the construction of your sampler it should be noted that your new sampler was fabricated and polished in a metal working factory. Please ensure that the sampler is thoroughly cleaned before the it is used for the first time.

If you have any questions regarding the use of this device please contact QAQC Lab