

MicroThief

General


The MicroMate is designed for low volume sampling of free flowing powders and granules.

Operation

1. Insert the sampler into the product, ensure that the tip is inside the sampler body.
2. At the required depth pull up the body to expose the tip. Powder will flow in around the tip.



3. Push down body of the sampler to trap the sample.
4. Withdraw sampler.
5. Pull up body to release the sample.

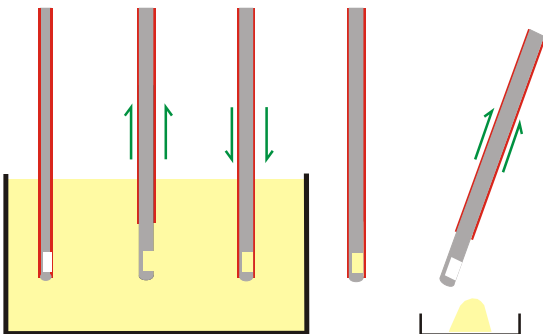


Cleaning

The sampler should be stripped down to its component parts and then cleaned according to the usual cleaning methods that are employed for items such as general dispensing ware, tablet punches etc.

Ensure that the all component parts are thoroughly dried after washing.


Hand Wash ✓	Washing Machine ✓	Ultra-sonic bath ✓	Autoclave 121°C ✓
-------------	-------------------	--------------------	-------------------

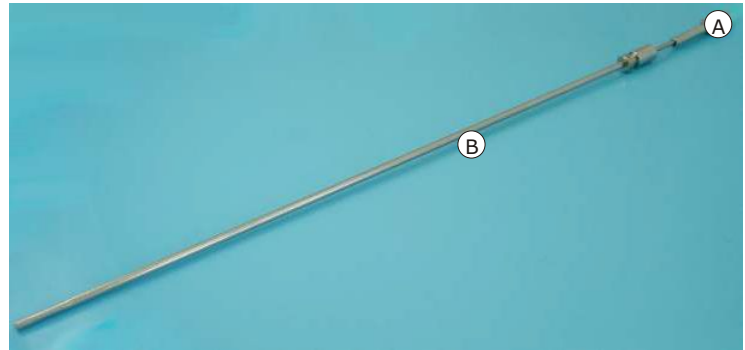

- (A) Removable handle
- (B) Outer tube

Adjusting The Sample Volume

- Loosen the retaining nut (E)
- Adjust the Sleeve
- Tighten the retaining nut (E)

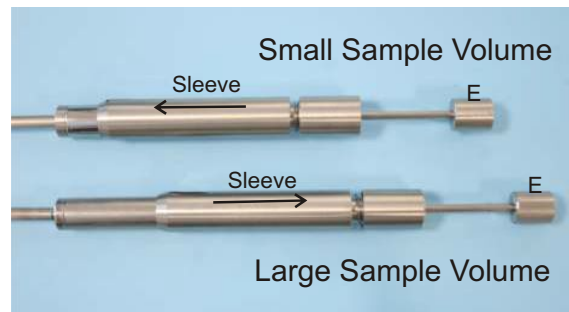

Note: You will have to test the sampler with your product to find the best position to achieve your required sample volume.

IMPORTANT: ENSURE SAMPLER IS CLEANED THOROUGHLY BEFORE INITIAL USE

Whilst every care has been made during the construction of your sampler it should be noted that your new sampler was fabricated and polished in a metal working factory. Ensure that the sampler is thoroughly cleaned according to your normal cleaning procedures and that any polishing waxes have been completely removed before the sampler is used for the first time.