

POWDER MAX


Large volume sleeve sampler

Designed for use with free flowing powders, granules and grains. The Powder Max enables large samples to be quickly and easily taken.

The Powder Max is made from 316 stainless steel.

Operation

1. Insert sampler into the product.
2. This action forces the sleeve upwards covering the sample hole and stopping powder from entering.
3. On pulling the sampler upwards, the sleeve moves down and the product can now enter the sampler.
4. Invert sampler to discharge the sample.

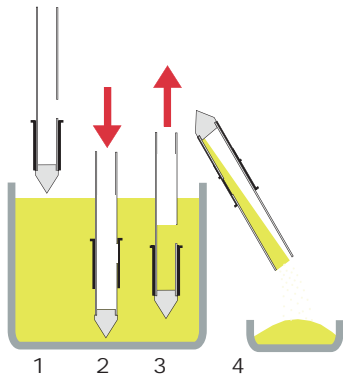

[QUOTE REQUEST](#) ✓

[BUY ONLINE](#) ✓

Stock No.	Length	Material	Diameter	Max Sample Vol.
675 1213A01	1000mm	316ss	40mm	150ml
Accessories	Cleaning brushes	See page 129		
	Storage Cases	See page 130		
	XRF Certification	See page 139		